



BIOPAC Systems, Inc.

Updated 01.04.07

42 Aero Camino
Goleta, CA 93117
Tel (805) 685-0066
Fax (805) 685-0067
www.biopac.com
info@biopac.com

Electrodermal Response (EDR) with an MP System

This Application Note will show you how to record galvanic skin response (GSR) and observe electrodermal response (EDR) changes with the galvanic skin response amplifier (GSR100C) or the telemetry (TEL100C) module of the MP System (with MP100 or MP150). With the MP System, you will measure both subtle and sudden changes in electrodermal activity, based on your protocols.

Electrodermal Response Basics

Electrodermal response (EDR) measurements show the activity of the eccrine sweat glands. Typically, one will place electrodes where the concentration of these glands is the highest: the fingertips. Responses are a function of the pre-secretory activity of eccrine sweat glands and the filling of the sudorific tubules. The combination of these sudorific elements serves to increase the conductivity of the skin when activated.

When one applies a feeble electric current (0.5V) between two electrodes (Ag-AgCl), the manifest electrical conductance varies in inverse proportion to the electric current flowing between the electrodes. The electrical conductance is a function of increasing eccrine activity. For instance, if a subject is presented a stimulus and the palms start to sweat, this response indicates a highly-stimulated state. The EDR of this subject will be higher than the subject's baseline. If another subject receives the same stimulus and the palms remain as "cool as a cucumber," the GSR reading remain unchanged with respect to the baseline. EDR undergoes relatively fast habituation (decrease of amplitude) in case the same stimulus is repeated over and over.

AcqKnowledge Software Features: EDR

- Automatically locate and analyze data
- Textual event markers
- On-line and off-line analysis
- Remote monitoring through the TEL100C
- On-line journal for note taking

- Automatic stimulation presentation (up to 16 on/off control channels)
- Export results to statistical programs for further analysis
- Determine maximum, minimum, and mean micromho (μmho) measurements

Applications: EDR

- Polygraph ("lie detector")
- Measure stress, arousal, emotional excitement
- Physiological biofeedback
- Measure absolute or relative response levels to different stimuli
- Relaxation Training

MP System Equipment: EDR

You can collect EDR data via the GSR100C or the TEL100C (the wireless model TEL100C-RF is not recommended for EDR applications).

- If using the **GSR100C**
 - Acquisition Unit: MP100 or MP150
 - Amplifier: GSR100C module
 - Transducer: TSD203 electrodermal response transducer set
 - Electrolyte: GEL101 skin conductance electrode paste or other electrolytic mixture
 - *Optional:* Module extension cable: MEC100C - 3 meters (extends lead length between GSR100C and TSD203)
- If using the **TEL100C**
 - Acquisition Unit: MP100 or MP150
 - Amplifier: TEL100C module
 - Transducer: SS3A electrodermal response transducer
 - Electrolyte: GEL101 skin conductance electrode paste or other electrolytic mixture

Hardware Setup

1. Set the **Gain** switch on the GSR100C.

To measure galvanic skin responses, you need to estimate the approximate units of the skin conductance (μmho) to be investigated.

- The unit of measurement for the GSR100C is the micromho

(μmho). The mho is the reciprocal of the unit of measurement for resistance, the ohm.

- Larger values indicate a higher level of conductivity; smaller values indicate less conductivity. For example:
 - If the subject is shown provocative slides, the response may be in the 0-50 μmho range.
 - If the subject is presented with a sudden 120dB buzzer in a quiet room, the response may be in the 0-200 μmho range.

Once you've determined which range will fit your experiment, choose the proper switch setting from the chart below. For example, for the 0-100 μmho range, set the module Gain switch to the 10 $\mu\text{mho/V}$ setting (on your module Gain switch, mho is an upside down ohm symbol).

Conductance Range		GSR100C Gain Switch
<i>DC</i>	<i>0.05 Hz</i>	
0-200 μmho	+/- 200 μmho	20 $\mu\text{mho/V}$
0-100 μmho	+/- 100 μmho	10 $\mu\text{mho/V}$
0-50 μmho	+/- 50 μmho	5 $\mu\text{mho/V}$
0-20 μmho	+/- 20 μmho	2 $\mu\text{mho/V}$

If using the [TEL100](#) Remote Monitoring System, for GSR measurements, the following gain settings correspond to mmhos. These settings can then be used to calibrate the signal using the rescaling feature in *AcqKnowledge* (see [AH-103](#) for more details on TEL100 setup).

TEL100 Gain	mmho/V
50	1,000
100	500
200	250

500	100
1,000	50
2,000	25
5,000	10
10,000	5
20,000	2.5
50,000	1

- Set the three **Filter** switches on the GSR100C to the appropriate filtering option for your protocol.

Low Pass (LP)	1 Hz LP	Sufficient for almost all EDR studies.
	10 Hz LP	Use to investigate higher frequency components of the EDR.
DC	DC	Gives direct (absolute) EDR readings from the subject.
High Pass (HP)	0.5 Hz HP	<u>Provides relative EDR recordings.</u> Removes low frequency signals. The effect of using the 0.5 Hz HP setting will be that your subject's baseline will return to nearly zero 1 second after a response change. The result provides an indicator of EDR changes as opposed to a specific EDR level.

0.05 Hz HP Provides relative EDR recordings. Removes very low frequency signals. The effect of using the 0.05 Hz HP setting will be that your subject's baseline will return to nearly zero 10 seconds after a response change. The result provides an indicator of EDR changes as opposed to a specific EDR level.

3. Set the **Channel** switch on the top of the GSR100C.
 - Select 1-16 to correspond with an available Analog Input channel.
4. Connect the **TS203** transducer to the GSR100C.
 - a. Plug the two colored electrode cables into the VIN+ and VIN- inputs.
 - Either blue lead can be connected to either VIN input.
 - b. Plug the black cable into the GND.

GROUNDING When using the GSR100C amplifier with other biopotential amplifiers attached to the same subject, do not attach the ground lead from the biopotential amplifier(s) to the subject. The subject is already appropriately referenced (grounded) to the system via the VIN- attachment to the GSR100C. If a biopotential ground is attached to the subject, then currents sourced from the GSR100C will be split to the biopotential amplifier ground lead, potentially resulting in measurement errors.

Optional: If using the module extension cable for up to three meters of extra distance between the subject and the MP acquisition unit, plug the **MEC100C** into the GSR100C inputs and then plug the TSD203 into the MEC100C inputs.

Software Setup

1. Open the **Change Scaling Parameters** dialog for the selected GSR100C channel (MP menu>Setup Channels>Scaling).
2. Enter **CAL1 Input Value** as 0.
3. Enter **CAL1 Scale Value** as 0.
4. Enter **CAL2 Input Value** as 1.
5. Enter **CAL2 Scale Value** to match the Gain setting on the GSR100C per the

chart below.

GSR100C Gain Switch	CAL2 Scale Value
20 $\mu\text{mho/V}$	20
10 $\mu\text{mho/V}$	10
5 $\mu\text{mho/V}$	5
2 $\mu\text{mho/V}$	2

- Enter the **Units** as μmho .
- Click **OK** to establish the settings and close out of the Scaling Parameters dialog.

For example... To use the 0-100 μmho range, set the switch on the GSR100C front panel to 10 $\mu\text{mho/V}$, set the channel on top of the module to the corresponding software channel, go to Scaling, and type in the following values for that channel:

Cal1: Input: 0, Scale 0; Cal2: Input 1, Scale 10; Units: μmho

The screenshot shows a dialog box titled "Change Scaling Parameters" for "A1, EDR (GSR100C)". It contains two rows of input fields: "Cal1" with "Input volts" set to 0 and "Scale value" set to 0; and "Cal2" with "Input volts" set to 1 and "Scale value" set to 10. The "Units" field is set to "mmho". There are "OK" and "Cancel" buttons on the right. At the bottom, there is an "Option" section with two unchecked checkboxes: "Calibrate ALL channels at the same time" and "Use mean value", and a "Settings..." button.


Scale settings for GSR100C @ 10 $\mu\text{mho/V}$ Gain

Optional Calibration

To verify the Gain setting of the GSR100C:

1. Calibrate *AcqKnowledge* for lower frequency response at DC:

- In the Scaling window, set the Input voltages so they map to the "DC" conductance ranges indicated by the sensitivity setting.
Cal 1: Input 0, Scale 0; Cal2: Input 1, Scale 5; Units: micromhos

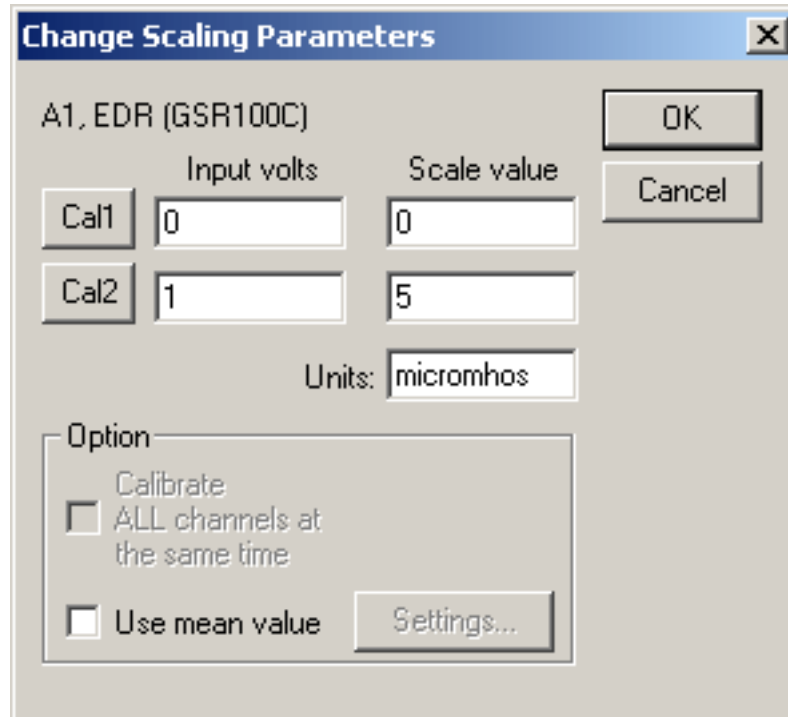

2. Place both lower frequency response (HP) filters on the GSR100C to **DC**.
3. Set the **Gain** switch on the GSR100C to 5 μ mho/V.
4. Perform measurement with electrodes disconnected.
 - *AcqKnowledge* should produce a reading of 0 μ mho. If slightly off, adjust the ZERO trim pot on the top front of the GSR100C to perfectly zero the reading.
5. Insulate a 100kohm resistor and place it from electrode pad to electrode pad (resistor must be insulated from fingers).
6. Perform measurement with electrode-resistor setup.
 - *AcqKnowledge* should produce a reading of 10 μ mho.

Subject Setup

The TSD203 transducer provides a small constant current to the site, and the measured resistance between the two electrodes constitutes the electrodermal response. When using the TSD203 to measure electrodermal response, you will need to determine the choice of electrolyte. For effective monitoring of local eccrine activity, use a higher impedance electrolyte with hyposaturated electrolyte concentrations of Cl⁻ (on the order of physiological levels). Many of our researchers have found electrolyte mixtures of 0.05M NaCl are optimal; BIOPAC GEL101 skin conductance electrode paste is a convenient option.

TSD203 Electrode Setup



GSR100C shown with MEC100C and TSD203

1. Apply GEL101 to the skin at the point of electrode contact and rub it in. The gel must have a chance to be absorbed and make good contact before recording begins.
2. Make sure the cavities of the TSD203 transducer are clean and fill them with GEL101 (or your own electrolyte mixture).
3. Attach the TSD203 electrode to the subject's fingertips as shown in the picture above, and then wrap the electrode around the fingers with the Velcro straps, applying mild tension.
4. Wait 5 minutes (minimum) before starting to record data.

[Return To Application Note Menu](#)